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The High-Performance Model with RES Technology

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**"It is a very robust mill that allows us to homogenize samples with a very uniform particle size according to the sieve used and very easy to use. Since it is a very robust equipment, as it is made to last, definitely I would recommend."**

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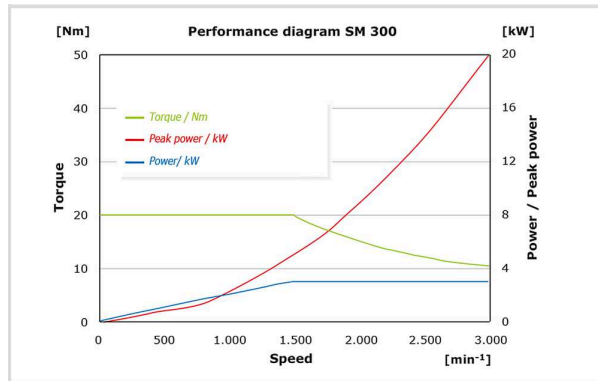
**OUR MOST POWERFUL CUTTING MILL FOR A WIDE RANGE OF APPLICATIONS**

This video shows the amazing versatility of the SM 300 cutting mill, with a wide range of accessories including different hoppers, collection vessels and rotors to facilitate the feeding and collection of sample materials of different shapes, sizes and quantities. The variable speed allows the perfect cutting parameters to be selected, and the powerful 3 kW motor with RES technology gives the mill maximum pulling power.

Watch the video to see how the SM 300 handles these 8 applications with flying colours:

- | Wood chips: Efficient size reduction of larger sample volumes
- | PC boards: Pulverising solid PCBs with the 6-disc rotor for precious metal analysis
- | Leather and table tennis balls: Comminution of textiles and rubber using the cyclone for rapid sample discharge from the grinding chamber
- | Secondary fuels: Efficient comminution of highly heterogeneous materials for rapid calorific analysis
- | Hay: Reproducible size reduction of fibrous, light and voluminous materials
- | Injection moulding sprues: Grinding at low speeds to minimise fines (dust) content
- | Toys: crushing of tough plastics by embrittling samples with liquid nitrogen

The SM 300 delivers perfect results even for the most demanding grinding tasks!

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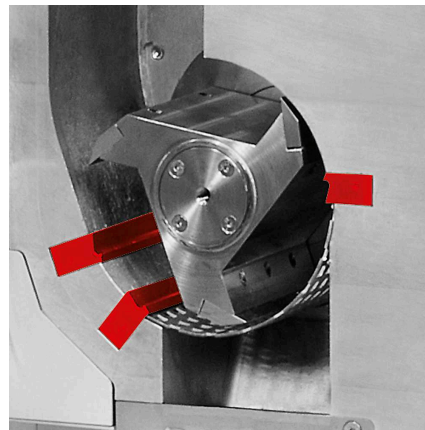
## TORREFACTION & BIO-COAL

Torrefaction is a thermochemical process carried out in the absence of oxygen to cause the partial decomposition of biomass. Water and volatiles are released from the biomass to leave a solid, dry material known as bio-coal. For some applications, bio coal can be used to replace traditional fossil fuels, thus helping to reduce CO<sub>2</sub> emissions.

In this video, both RETSCH cutting mill SM 300 and the CARBOLITE rotating tube furnace TSR have been used as a part of a torrefaction process to create bio-coal fuel.

The Verder Scientific companies provide solutions to support your process every step of the way, from sample preparation and analysis, to heat treatment, and more.

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**Cutting Mill SM 300 316L complete with collecting receptacle, 5 litres (316L) and base frame with wheels (please order rotor, hopper and bottom sieve separately)**

20.729.0010  SM 300, 220–230 V, 50/60 Hz, 속도 변환 용 주파수 변환기 포함, housing and double acting cutting bars of steel 316L, foodGrade

20.729.0011  SM 300, 220–230 V, 50/60 Hz, 속도 변환 용 주파수 변환기 포함, housing of steel 316L, foodGrade, double acting cutting bars of FDA certified steel, foodGrade

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#### ROTOR, STEEL 316L, FOODGRADE

22.608.0037 Parallel section rotor, cutting plates of steel 316L, foodGrade

22.608.0038 Parallel section rotor, cutting plates of FDA certified steel, foodGrade

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22.020.0009 Cyclone unit (316L) incl. collecting receptacle, 30 litres, steel 316L and carriage  
for collecting receptacle (to be used with industrial vacuum cleaner)

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ACCESSORIES FOR PARALLEL SECTION ROTOR

22.151.0009 Cutting plates, steel 316L, for parallel section rotor, 1 set (3 pieces)

22.151.0008 Cutting plates, FDA certified steel, for parallel section rotor, 1 set (3 pieces)

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22.152.0011 Double acting cutting bars, steel 316L, 1 set (3 pieces) for SM 300 316L,  
foodGrade

22.152.0010 Double acting cutting bars, FDA certified steel, 1 set (3 pieces) for SM 300 316L,  
foodGrade

## ADDITIONAL ITEMS SM 300 316L, FOODGRADE

99.200.0014

SM 300 IQ / OQ 문서