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**The RETSCH AS 200 tap is suitable for dry sieving with test sieves of 200 mm or 8" diameter. The combination of horizontal, circular sieving motions with vertical taps reproduces the principle of hand sieving. The uniform mechanical action ensures reliable and reproducible measurement results.**

This special type of sieving motion used by the AS 200 tap is specified in various standards for particle size analysis of materials such as activated carbon, diamonds, spices, metal powders, abrasives or cement. Operating the AS 200 tap is exceptionally easy and safe. The integrated clamping device allows for sieve stacks with up to 7 or 13 fractions, depending on the height of the sieve frame. The sieving time is set from 1 to 99 minutes via a digital display. The number of rotations and taps is fixed; the tapping motion can be deactivated, if required. A safety switch and an anti-trap protection provide maximum safety. Thanks to an integrated interface, the AS 200 tap can be controlled with the evaluation software EasySieve<sup>®</sup>.

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- | Sieving with circular motion and vertical taps according to standards
- | Measuring range 20 μm to 25 mm
- | For 200 mm / 8" sieves
- | Sieve stack up to 350 mm
- | Digital time setting
- | Integrated interface
- | Suitable for dry sieving

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RETSCH sieve shakers can be equipped with a variety of accessories to meet a wide range of application requirements.

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**RETSCH TEST SIEVES AND ACCESSORIES - ENGINEERED FOR SUPERIOR PERFORMANCE**

## SOFTWARE DE EVALUARE EASYSIEVE / EASYSIEVE CFR

EasySieve, the software for particle size analyses, exceeds manual evaluation in many aspects. The software is able to automatically control the necessary measurement and weighing procedures – from the registration of the weight of the sieve up to the evaluation of the data. It is simple and convenient to use and is also available in an FDA 21 CFR Part 11-conform version.

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## PROBE TIPICE

RETSCH's Analytical Sieve Shaker AS 200 tap is ideally suited for separation, fractioning and particle size determination of abrasives, charcoal, cement, coal, diamonds and spices.

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[www.retsch.ro/as200tap](http://www.retsch.ro/as200tap)

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### TAP SIEVE SHAKER AS 200 TAP

**Tap Sieve Shaker AS 200 tap for test sieves 200 mm / 203 mm (8") Ø, incl. sieve lid and EasySieve evaluation software (please order test sieves and collecting pan separately)**

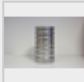
30.025.1001  AS 200 230 V, 50 Hz incl. sound-enclosure cabinet, conforms with CE standards  
tap

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### SIEVE STACKS AS 200 TAP

60.131.000998  Sieve stack consisting of 6 test sieves (ISO 3310-1), 200 mm Ø, 50 mm height (63 µm, 125 µm, 250 µm, 500 µm, 1 mm, 2 mm) and collecting pan

60.150.000998  Sieve stack consisting of 6 test sieves (ASTM E11), 203 mm (8") Ø, 50 mm (2") height (230 mesh, 120 mesh, 60 mesh, 35 mesh, 18 mesh, 10 mesh) and collecting pan

### ADDITIONAL ITEMS AS 200 TAP

22.783.0001  Ball-pan hardness test kit acc. to ASTM D3802-79

99.200.0021 IQ/OQ Documentation for AS 200 tap

FOR TEST SIEVES 200 MM Ø, HEIGHT 50 MM

69.720.0050  Collecting pan oțel inoxidabil 200 mm Ø height 50 mm

FOR TEST SIEVES 200 MM Ø, HEIGHT 25 MM

69.720.0025  Collecting pan, oțel inoxidabil, 200 mm Ø, height 25 mm

FOR TEST SIEVES 203 MM Ø / 8" Ø, HEIGHT 2"

69.720.3050

Collecting pan, oțel inoxidabil, 8" Ø, height 2"

FOR TEST SIEVES 203 MM Ø / 8" Ø, HEIGHT 1"

69.720.3025

Collecting pan, oțel inoxidabil, 8" Ø, height 1"